



149 Entin Road  
Clifton, New Jersey 07014  
973.594.1100  
973.594.1500 - Fax

## HSLT Series Science Lab Table Leg Assembly Instructions

CLIFTON NJ 07014  
ONTARIO CA 91764  
MEMPHIS TN 38106

The HSLT Series has been designed with the capability to be used with both our standard non- book compartment style or with book compartments (purchased separately).

To assemble the table using the standard non-book compartment style, follow the below instructions.

### Tools Required

WRENCH



**Step 1:** Place the tabletop upside-down on the floor with apron side up.

*Note: Floor should be clear of any objects or dirt in order not to damage the tabletop.*

**Step 2:** Remove plastic covers on bolts and place each leg in position with leg bolts going through the corner braces.

**Step 3:** Place a washer, lock washer and nut onto each bolt and tighten well using the provided wrench.

**Step 4:** Slide a leg boot onto each leg assuring that the wider side is facing the bottom of the leg.

**Step 5:** Screw a glide into the bottom of each leg.

**Step 6:** Turn the table over with the tabletop facing up. Place the table in desired location. Adjust the glides to assure that table is stable.

**Step 7:** Slide down protecting boots to cover the adjusting glides.

*Note: Extra apron marked BC is for use with book compartments only.*

---

Dear Customer,

Thank you for your purchase. Although extreme care goes into manufacturing & packaging of this product, mistakes are possible. If you encounter any missing parts or difficulty in assembly. etc. Please call our customer service **TOLL FREE at 1-800-261-4112** for courteous & immediate resolution to your problem.

**CAUTION:** Double check all screws to be tightened securely before use. Check all screws periodically.