

Technical Information Sheet

Product Number 9792R

3M[™] Aluminum Tape for Diagnostic and Bioanalytical Constructions

Effective: May 2010

Supersedes all previous specifications.

Product Description:

3M[™] Aluminum Diagnostic Tape 9792R consists of a "dead soft" aluminum foil for high conformability coated on one side with a pressure sensitive acrylate adhesive. The tape is supplied on a brown silicone coated paper release liner. Liner wound "IN".

	plications: Bioanalytical devices		Microplate Seals
Fe	atures and Benefits:		
	Highly Conformable		Low level of extractables in a variety of solvents
	Pierceable		Adheres well to a wide range of substrate
	Suitable for use between -70°C and 100°C (-94°	F ar	nd 212°F)
<u>ر</u> م	mnosition:		

Composition:

Backing	Aluminum Foil	1.4 mils (0.04 mm) typical
Adhesive	Pressure sensitive acrylate, neutral	1.1 mils (0.03 mm) typical
Release Liner	61# Brown Kraft Paper, one-side silicone coated	4.6 mils (0.12 mm) typical

Typical Product Properties¹

Typical Froduct Froperties.	
Adhesion to Steel ²	35 oz/in [1.1 kg/25 mm] [9.7 N/25 mm]
Liner Release ²	10 g/in [10 gms/25 mm] [0.1 N/25 mm]
Solvent Extractables (Reference: GC/MS) 80% Ethanol/20% Water:	None detected above background (typical) 5 μg/ml hydrocarbon acrylate esters (typical) 630 μg/ml antioxidant (typical)

¹ 3M test methods are available upon request.

Product Dimensions: Product supplied on 3-inch (76 mm) white plastic cores.

Roll Length, Maximum	500 yd (457 m) ± 2%
Roll Width, Maximum	42 in (106.7 cm) ± 1/16 in (1.6 mm)

Splice Frequency: Intended for continuous processing.

Please contact your 3M Medical	Roll Length:	Maximum # of Splices:
Specialties Account Representative for	1 to 100 yd (0.9 to 91 m)	2
	101 to 200 yd (92 to 183 m)	3
information on	201 to 400 yd (184 to 366 m)	5
splice-free rolls.	401 to 500 yd (367 to 457 m)	7

Recommended Storage Conditions & Shelf Life: Product as supplied in original packaging will maintain stated test properties for a period of two years from date stamped on shipping container when stored at temperatures of 50-80°F (10-27°C) and relative humidity of 40-60 percent.

Please see last page for important product, safety and warranty information.

² Lot specific results may be included on a Certificate of Analysis upon request at time of order.

Product and Safety Information: User is solely responsible for determining the suitability of 3M samples and products for the intended use including any necessary safety or toxicity assessment. 3M will provide Material Safety Data Sheets and results of toxicity testing upon request. In every case before using any product in full scale production users should conduct their own tests to determine to their own satisfaction whether the product is of acceptable quality and is suitable for their particular purposes under their own operating conditions.

Notice: Nothing contained herein shall be construed to imply the nonexistence of any relevant patents or to constitute a permission, inducement or recommendations to practice any invention covered by any patent, without authority from the owners of this patent.

All statements, technical information and recommendations herein are based on tests 3M believes to be reliable, but the accuracy or completeness thereof is not guaranteed. 3M warrants only that products will meet 3M's specifications at the time of shipment to the customer. 3M does not offer any other warranty and does not warrant the performance, safety or such other characteristics of Products in combination with other materials. 3M specifically DOES NOT warrant Products for any intended or unintended uses (whether or not foreseeable); for compatibility or suitability with other components or compatibility with any methods of manufacture or conversion. The foregoing warranty is made in lieu of all other warranties, expressed or implied, including the implied warranties of merchantability, fitness for a particular purpose and freedom from non-infringement.

Limitation of Remedies: If products are proven not to meet 3M's specifications, the sole and exclusive remedy available and 3M's only obligation shall be, at 3M's option, to replace such quantity of Products which are proven out of specification or to refund the purchase price paid for Products.

Limitations of Liabilities: The remedies provided herein are exclusive remedies against 3M for any alleged or actual nonconformance to specifications or defect or other failure in products or for 3M's performance of its supply obligations. Under no circumstances is 3M liable for any direct, indirect, incidental, special or consequential damages (including lost profits) in any way related to the product under any theory of law including, but not limited to, negligence and strict liability.

Ordering Information

Call us at Customer Service to place an order: 800-742-1994 (U.S.). Visit our website: www.3M.com/medicalspecialties for product and services information, news about conferences we will be attending, new product highlights or to make a direct inquiry.

To have a sales representative contact you, to request samples or clinical and safety summaries, please contact us at 3M HELPLINE 800-228-3957 (U.S.) or for international inquiries, please contact your local country representative. Our 3M Medical Specialties subsidiary contacts are listed below.

International Locations

Australia 3M Health Care Tel: 1 300 363 878 Fax: 1 800 060 888

Belgium 3M Belgium N.V./S.A. Tel: 32-2-722-5111 Fax: 32-2-720-0225

Brazil 3M do Brasil Ltda. Tel: 55-19-3838-6661 Fax: 55-19-3838-7449

Canada
3M Canada
Tel: 800 364-3577
Fax: 800 341-4630

Chile 3M Chile S.A. Tel: (56-2) 4103000 Fax: (56-2) 4103400

China 3M China Ltd. Tel: 86-21-6275-3535 Fax: 86-21-5208-2205

Colombia 3M Colombia S.A. Tel: 57-1 4101666

Fax: 57-1 4161677 Czech Republic

3M Cesko s.r.o. Tel: +420-261 380 111 Fax: +420-261 380 110

Denmark 3Ma/sTel: 45-43-48-0100

Fax: 45-43-96-8596

Argentina **Dominican Republic** 3M Argentina S.A.C.I.F.I.A. 3M Dominicana Tel.: 54-11-4339-2400 Tel: (809) 530-6560 X499 Fax: 54-11-4339-2640 Fax: (809) 530-2960

Finland Suomen 3M Oy Tel.: 358-9-52-521 Fax: 358-9-512-29-44

France	
Laboratoires 3M San	té
Tel: 33-1-3031-8376	,
Fax: 33-1-3031-8378	

Hong Kong
3M Hong Kong Ltd.
Tel: (852) 2806 6111
Fax: (852) 2234 6044

Hungary
3M Hungária Kft.
Tel: (36-1) 270-7777
Fax: (36-1) 320-0951

India
3M India Ltd.
Tel: +91-80-2223 1414
Fax: +91-80-22231431

Indonesia
PT 3M Indonesia
Tel: (62-21) 520 3401
Fax: (62-21) 520 3106

Israel	
3M Isra	ael Ltd.
Tel: 97	2-9-9615000
Fax: 97	2-9-9615050

Italy
3M Italia SPA
Tel: +39 2 7035-2406
Fax: +39 2 7035-2484

Janan 3M Health Care Ltd. Tel: 81-3-3709-8289 Fax: 81-3-3709-8754

Kazakhstan
3M Representation Office
Tel: +7 3272 509 944
Fax: +7 3272 509 573

Korea
3M Korea Ltd.
Tel: 82-2-3771-4286
Fax: 82-2-786-2825

Malaysia
3M Malaysia Sdn. Bhd
Tel: 6-03-7806 2888
Fax: 6-03-7806 2902

Mexico
3M México, S.A. de C.V
Tel: +52-55-5270-0400
Fax: +52-55-5270-0433

Middle East
3M Gulf Ltd.
Tel: 971-4-3670-777
Fax: 971-4-3670-699

The Netherlands
3M Nederland B.V.
Tel: 31-715-450-450
Fax: 31-715-450-212

New Zealand
3M New Zealand
Tel: 64-9-477 4040
Fax: 64-9-477 6692

Peru
3M Peru S.A.
Tel: 511-224-2728
Fax: 511-224-3171

Philippings

Poland 3M Poland Sp.z o.o. Tel: (48 22) 739 60 00 Fax: (48 22) 739 60 01

Thailand

Taiwan

Turkey

Ukraine

3M Ukraine

Venezuela

Viet Nam

3M Turkey

3M Thailand Ltd.

3M Taiwan Ltd.

Fax: (66)2 261 7535

Tel: 886-2-2704-9011

Fax: 886-2-2706-0355

Tel: 9 (0212) 3507777

Fax: 9 (0212) 2821742

Tel: +380 44 490 57 77

Fax: +380 44 490 57 75 **United Kingdom & Ireland** 3M Health Care Ltd. Tel: 44-1509-611-611

Fax: 44-1509-237-288

Tel: 58-241-8391911

Fax: 58-241-8391928

3M Viet Nam Ltd.

Tel: 84 8 640986

Fax: 84 8 641024

3M Manufacturera Venezuela

Tel: (66)2 260 8577 Ext.175

Portugal
3M Portugal
Tel: 351-21-3134500
Fax: 351-21-3134680

Russia	
3M Russia	
Tel: 7 (495) 784	-74-74
Fax: 7 (495) 784	1-74-75

Singapore
3M Technologies (S) Pte. Ltd
Tel: 65-6450-8866
Fax: 65-6455-2130


Sri Lanka
3M Lanka (Pvt.) Ltd.
Tel: (94) 11 2785701
Fax: (94) 11 4410083-4

Spain
3M España, S.A.
Tel: 34-91-321-6000
Fax: 34-91-321-6002

Sweden
3M Svenska AB
Tel: +46-8-92 21 00
Fax: +46-8-92 24 22

Switzerland & Eastern Europe

3M (Schv	vei	z) A	G	
Tel:	+41	44	724	92	44
г	. 41	4.4	724	0.4	00

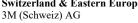

Fax: +41 44 724 94 80



3M Medical Specialties

3M Center, Building 275-5W-05 St. Paul, MN 55144-1000 USA Tel.: 800 228-3957